**Product number: K8-7035** 

Product name: Seta-750-mono-NHS

Date: 05/24/11 Page 1

#### **General Data**

Molecular Mass: 1268.58

1010.10 (protonated form)

**Solubility:** Water, Alcohol, DMF, DMSO **Insoluble:** Acetone, Chloroform, Toluene

Storage: Store in absence of light, desiccate and refrigerate

# **Description**

Seta-750-mono-NHS (K8-7035) is a hydrophilic, amine-reactive label containing one NHS-ester group with almost identical absorption and emission as Cy7 or Alexa 750 and can therefore be used with these filter sets. It combines high photostability and brightness. Its extinction coefficient is  $235,000 \, \mathrm{M}^{-1} \mathrm{cm}^{-1}$ .

## **Applications**

- · Covalent labeling of proteins, amino-modified DNA and amino-modified oligonucleotides
- Fluorescence intensity and fluorescence polarization-based applications
- Resonance Energy Transfer (RET)
- Flow Cytometry
- Immunofluorescence
- · Gene Expression
- Homogeneous Assays
- Microarrays

#### **Advantages**

- Perfectly suited for excitation with the 695 nm or 740 nm diode lasers
- Sensitive; high extinction coefficients and high quantum yields
- pH-insensitive between pH 3 and pH 10
- Good aqueous solubility; this label does not alter the solubility of bioconjugates
- High photostability; e.g. compared to fluorescein or Cy7
- Low molecular weight Seta-750 does not add substantial mass to the conjugates
- · Ideal for non-radioactive labeling of proteins, amino-modified oligonucleotides and amino-modified lipids

## **Spectral Data**


Solvent System: phosphate buffer pH 7.4

Sample	Dye-to-protein Ratio	Absorption max. [nm]	Extinction Coefficient [M <sup>-1</sup> ·cm <sup>-1</sup> ]	Fluorescence* max. [nm]	Quantum Yield [%]
Free dye	-	751	235,000	779	13
IgG conjugate 1	1.0	756		782	14
IgG conjugate 2	2.0	755		782	12
IgG conjugate 3	4.0	755		782	10


<sup>\*</sup> Excitation at 700 nm

**Product number: K8-7035** 


Product name: Seta-750-mono-NHS

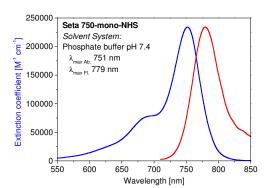


Absorption and emission spectrum of **Seta 750** in phosphate buffer (pH 7.4)

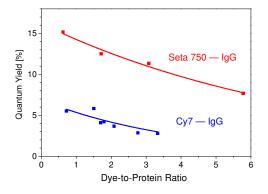

Quantum yield vs. dye-to-protein ratio (D/P) of **Seta 750 — IgG** conjugates in phosphate buffer (pH 7.4)

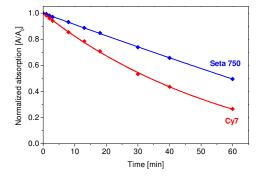


Decrease of the long-wavelength absorption of **Seta-750** compared to **Cy7** upon irradiation with a Metalhalogen lamp Philips HPI-T Plus 400W/645

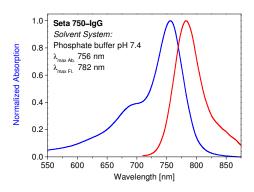


Date: 05/24/11 Page 2

Absorption and emission spectrum of a **Seta 750 — IgG conjugate** in phosphate buffer (pH 7.4, Dye-to-protein ratio ~0.6)

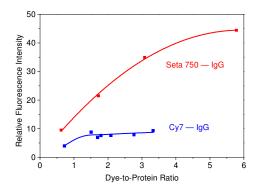

Relative fluorescence intensity (Q.Y. × D/P) vs. dye-to-protein ratio (D/P) of **Seta 750 — IgG conjugates** in phosphate buffer (pH 7.4)

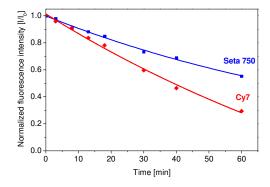

Decrease of fluorescence intensity of **Seta-750** compared to **Cy7** upon irradiation with a Metalhalogen lamp Philips HPI-T Plus 400W/645